

A complete, fully programmable blot processing station



The Freedom Rocker™ BlotBot® Automated Blot Processor

KEY BENEFITS

► IMPROVE CONSISTENCY

Microprocessor controls ensure precise timing and dispensing. Get the same great results every time you use it.

► CONSERVE ANTIBODY

Only 3.5 mL of antibodies needed for each miniblots. 3 mL can be recovered.

► USE EXISTING PROTOCOLS

No need to change the way you process your blots. Programming is easy with a color graphic display.

“I’m one of the first users of the [previous] Freedom Rocker at Amgen in Cambridge, MA. The machine gets LOTS of use from all of us here.”

DR. BARBARA GRUBINSKA, AMGEN

Stop babysitting your Westerns! Incubate and wash your samples unattended over night or during a busy day with the next generation of Freedom Rocker.

Save Time and Money

This fully programmable blot processing station automates the tedious incubation and washing steps to process blots and gels, saving you time and money through increased productivity. You can be confident that all steps will be performed with consistency. All trays receive the same reagent or wash buffer at the same time in equal volumes.

Simply dispense individual reagents (as little as 3.5 mL) into each tray and recover them separately. The Freedom Rocker is perfect for conserving precious primary antibodies.

User Friendly Interface

The control panel has a large graphic display and an intuitive layout, making it very easy to program your own protocols, which can be very complex with many steps. Programs can include different rocking speeds and times for each step.

Proven Track Record

Freedom Rockers are built to last. In fact, many of the original Freedom Rockers from 2005 are still in use today.

Processing a Western Blot



The Freedom Rocker frees up your time, automating dozens of liquid handling steps and ensuring consistency.

Features

- Automated delivery and removal of buffers/reagents
- Dispense and recover different antibodies or reagents in each tray
- Specially designed trays require as little as 3.5 mL of reagent per blot
- Reagent recovery for added cost savings
- The Freedom Rocker™ can call any lab bench home. Just plug into an electrical outlet
- Quiet — no air connections or vacuum pumps required
- Drip proof quick-release connections
- Patent pending technology
- 2 year warranty
- 30-day money back guarantee

“The Freedom Rocker is perfect for silver staining of gels for mass spectrometry. It has made an unbelievable impact on data generation for me in cancer protein screening. “

DR. STEVE GRIFFITHS, ATLANTIC CANCER RESEARCH INSTITUTE

“The blot washer you sent us is perfect! I love it!”

RENEE CLARK, CANADIAN FOOD INSPECTION AGENCY, LETHBRIDGE AB



For More Information, Contact: